


Chlorinated compounds on ZB-MR2**Column:** Zebtron[™] ZB-MultiResidue[™]-2, GC Cap. Column 30 m x 0.25 mm x 0.20 μ m, Ea**Phase:** Proprietary Pesticides Phase**Dimensions:** 30 meters x 0.25 mm x 0.2 μ m**Order No:** 7HG-G017-10**Oven Profile:** 40 °C for 2 min to 260 °C @ 15 °C/min to 280 °C @ 8 °C/min for 3 min to 320 °C @ 15 °C/min for 5 min.**Carrier Gas:** Constant Flow Helium, 1.5 mL/min**Injection:** Splitless :1 1 μ L @ 240°C**Detection:** Mass Selective (MSD) (320°C)**Analyst Note:** Analytes are 50ppm in Dichloromethane**Products used in this application:**

Chlorinated compounds on ZB-MR2**ANALYTES:**

- 1** 1,2-Dichlorobenzene
- 2** bis(2-Chloroethyl)ether
- 3** 1,4-Dichlorobenzene
- 4** 1,3-Dichlorobenzene
- 5** bis(2-Chloroisopropyl)ether
- 6** Hexachloroethane
- 7** bis(2-Chloroethoxy)methane
- 8** Hexachlorobutadiene
- 9** Hexachloropropene
- 10** 1,2,4-Trichlorobenzene
- 11** Hexachlorocyclopentadiene
- 12** 1,2,4,5-Tetrachlorobenzene
- 13** 2-Chloronaphthalene
- 14** Pentachlorobenzene
- 15** 1,3-Dinitrobenzene
- 16** 4-Chlorophenylphenylether
- 17** 4-Bromophenyl phenyl ether
- 18** Hexachlorobenzene
- 19** Pentachloronitrobenzene
- 20** Isodrin
- 21** Aramite
- 22** Aramite
- 23** Chlorobenzilate
- 24** Kepone

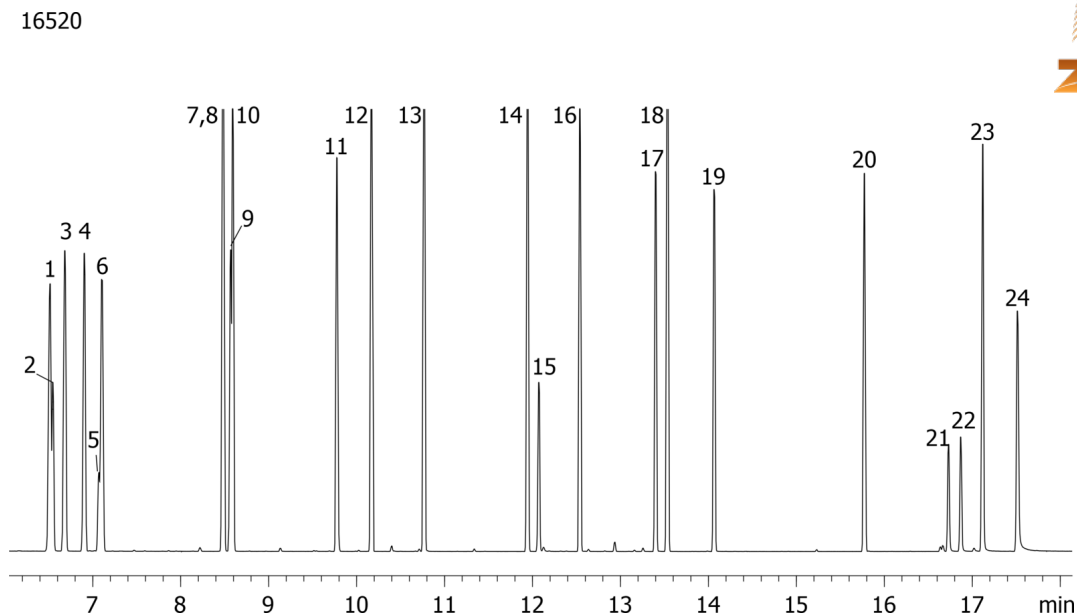